CertoClav Vacuum Pro Series LABORATORY AUTOCLAVE





CertoClav Vacuum Pro Series SUITABLE FOR ANY APPLICATION





CertoClav Vacuum Pro Series STAND-ALONE OR SYSTEM INTEGRATED



AUTOCIAVE ORDER CONFIGURATIONS

	12 liters	18 liters	22 liters	tanks (Stand alone)	connection. Including R/O- System for automatic refill
✓				✓	
✓					✓
	✓			√	
	√				√
		√		√	
		√			✓
			✓	√	
			✓		√
					alone) / / / / /







Reverse osmosis system Automatic water supply for the autoclave. Requires tap water connection. Power: 110-220V, 50-60Hz. Order number: 8500499




Label printer Print expiration date and protocol information on self-adhesive label and mark your goods. Order number: 8500502

Service set

Includes door sealing, air filter, safety valve, quick-connect drain hose Order number: 8500495

ACCESSORIES



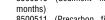

Order numbers:

Filters for reverse osmosis system

8500510 (Sediment filter, replace every 6 8500511 (Precarbon filter, replace every 6 months) 8500512 (Membrane filter, replace every 30 months)


Tray for bottles

Order numbers: 8I: 8500478A 12I: 8500479A 18I: 8500480A 22I: 8500481A

Barcode scanner Track the complete sterilization history of an item.

Order number: 8500503


External printer Print protocols with a printer Order number: 1300187

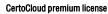

- Enable SMS notifications
- Storage for 100.000 protocols •
- Premium Remote support
- Real-time-data online
- Order number: 8500212

CertoClav Vacuum Pro Series INTUITIVE USER INTERFACE

TOUCH SCREEN

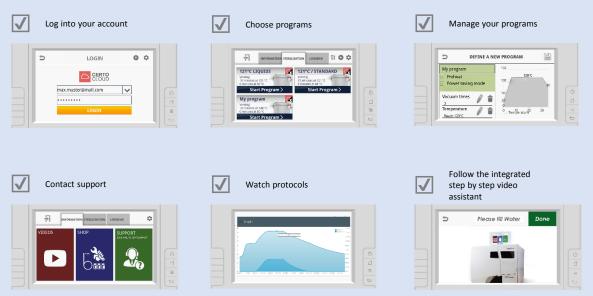




FREE AUTOCLAVE MANAGEMENT SOFTWARE

Access real-time data



List of Protocols				
Cycle 个	Program Name	Start		
76	cloud generated	7/22/ PM		
75	121°C SOLID	7/22/ PM		
74	Neues Programm	7/22/ PM		
74	121°C SOLID	7/2. PM		



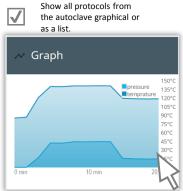




	with a few clicks.	,0013
Pro	otocol Verification	
₿	Protocol ID *	
-	Device Key *	0 / 24
		0/17
	Security Code *	01640
	VERIFY	

CertoClav Vacuum Pro Series STERILIZE SOLID MATERIALS AND LIQUIDS

STERILIZATION OF SOLID MATERIALS

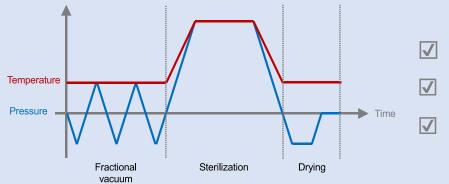


All sterilization programs for solids follow a typical Class-B cycle.

Precise combination of vacuum, pressure and temperature for optimal sterilization of solids.

The fractionated vacuum presses the saturated water vapor through the sterilization package and into each microscopic gap.

STERILIZATION OF LIQUIDS WITH MEDIA SENSOR

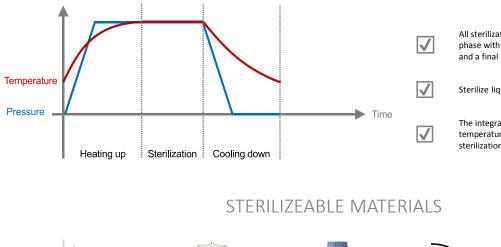

All sterilization programs for liquids follow a heating phase with venting followed by a sterilization phase and a final cooling phase..


Sterilize liquids at 121°C for 20 minutes.


The integrated media sensor measures the real temperature of liquids and ensures an optimum sterilization result.



Instruments

Textiles

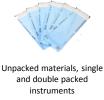

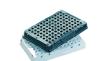

Plastics, rubber materials



Liquids in bottles







Liquids with media temperature sensor



Hollow materials



Foodstuffs



Filters, waste

CertoClav Vacuum Pro Series CUSTOM PROGRAM AND AUTOMATED MATERIAL TESTS

	Vacuum times	Temperature [°C]	Holding time [min]	Drying time [min]	Total time [min]
Solid	1	121	20	3	32
Solid quick	1	134	4	3	20
Porous	3	121	20	7	50
Porous quick	3	134	4	7	27
Hollow	3	121	20	10	47
Hollow quick	3	134	4	10	34
Liquid	0	121	20	0	45 to 120
Prion	3	134	20	10	50
B-D test	3	134	4	7	31
Vacuum test	1	-	15	-	15
Clean process	-	-	-	-	29
User-Defined	1 to 10	105 to 134	1 to 60	1 to 40	

PROGRAMS


MATERIAL TESTS





CertoClav Vacuum Pro Series STERILIZE SOLID MATERIALS AND LIQUIDS

LOADING CAPACITY

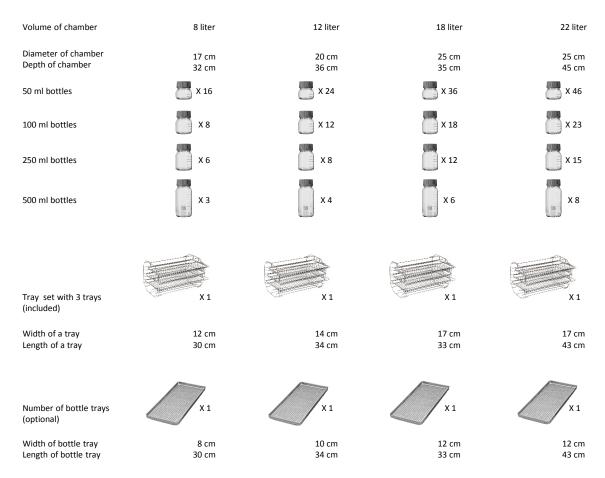



CertoClav Vacuum Pro Series THE ALL-IN-ONE AUTOCLAVE

TECHNICAL DATA

Chamber volume	8 liter	12 liter	18 liter	22 liter
Product category	Laboratory horizontal table-top autoclave	Laboratory horizontal table-top autoclave	Laboratory horizontal table-top autoclave	Laboratory horizontal table-top autoclave
Order number for autoclave with integrated water tanks (stand-alone)	8500486	8500487	8500488	8500489
Order number for autoclave with fixed water connection. Including RO-System for automatic refill	8500486W	8500487W	8500488W	8500489W
Vacuuming	√	√	\checkmark	√
Drying	√	√	√	√
Liquid media sensor	√	√	√	√
User-defined program	√	√	√	√
Printer port	√	√	√	√
Ethernet	√	√	√	√
WiFi	√	√	√	√
CertoCloud Software	√	√	√	√
Product size	56 x 44 x 36 cm	61 x 50 x 43 cm	61 x 50 x 43 cm	61 x 50 x 43 cm
Diameter of chamber	17 cm	20 cm	25 cm	25 cm
Depth of chamber	32 cm	36 cm	35 cm	45 cm
Net weight	35 kg	51 kg	53 kg	55 kg
Gross weight	42 kg	59 kg	61 kg	63 kg
Carton size	65 x 45 x 42 cm	73 x 57 x 51 cm	73 x 57 x 51 cm	73 x 57 x 51 cm
Scope of delivery	1 autoclave, 3 trays, 1 tray holder, 1 tray handle, 1 power cable, 1 quick-connect drain hoses, 1 CertoCloud license	1 autoclave, 3 trays, 1 tray holder, 1 tray handle, 1 power cable, 1 quick-connect drain hoses, 1 CertoCloud license	1 autoclave, 5 trays, 1 tray holder, 1 tray handle, 1 power cable, 2 quick-connect drain hoses, 1 CertoCloud license	1 autoclave, 3 trays, 1 tray holder, 1 tray handle, 1 power cable, 1 quick-connect drain hoses, 1 CertoCloud license
Sterilizable materials	Solids, hollow materials, porous materials, liquids, prions, textiles, plastics	Solids, hollow materials, porous materials, liquids, prions, textiles, plastics	Solids, hollow materials, porous materials, liquids, prions, textiles, plastics	Solids, hollow materials, porous materials, liquids, prions, textiles, plastics
Locking mechanism	Electric door lock	Electric door lock	Electric door lock	Electric door lock
Cooling system	Air cooling	Air cooling	Air cooling	Air cooling
Screen	7" Touch screen	7" Touch screen	7" Touch screen	7" Touch screen
Chamber material	Stainless steel 304	Stainless steel 304	Stainless steel 304	Stainless steel 304
Protocol export	USB, SD-card, cloud	USB, SD-card, Cloud	USB, SD-card, cloud	USB, SD-card, cloud
Languages	English, German	English, German	English, German	English, German
Sterilization temperature	121°C to 134°C	121°C to 134°C	121°C to 134°C	121°C to 134°C
Maximum working pressure	230 kPa	230 kPa	230 kPa	230 kPa
Power supply	220V, 50/60Hz, 2400W	220V, 50/60Hz, 2400W	220V, 50/60Hz, 2400W	220V, 50/60Hz, 2400W
Certification	CE	CE	CE	CE
Warranty	2 years	2 years	2 years	2 years



Certoclav Sterilizer GmbH Georg Grinninger-Str. 37 A-4050 Traun



Phone: +437229689 Email: support@certoclav.com Website: www.certoclav.com



Orders, dealer verification Custom manufacture Operating instructions and operation videos Technical support

```
Catalog 2017
```